

Exploring the role and measurement of particle size and shape in metal additive manufacturing

Introduction

In recent years, metal additive manufacturing (AM) has transitioned successfully from a prototyping tool to a still new, but established and economically viable, choice for component production. Indeed, as the aerospace, energy, automotive, medical, and tooling industries have embraced this technology [1], annual sales of metal AM machines have risen from less than 200 in 2012 to almost 2,300 in 2019. Alongside this trend, the use of AM in manufacturing is also driving an increase in the proportion of the market dedicated to metal materials. This segment is expected to account for a quarter of the whole manufacturing market by 2023 [2].

Powder bed AM processes such as selective laser melting (SLM), electron beam melting (EBM), and binder jetting offer certain advantages over alternative powder metallurgy methods, including design flexibility and potential for high material-use efficiency. These processes are particularly suitable for producing small to medium volumes of relatively small components, as well as enabling the creation of new, complex parts that were previously unachievable. As these technologies are adapted to enable larger components and higher throughputs, the development of AM machines is an increasingly important focus area. However, there is now also an equal emphasis on the properties of the powders used [3].

What makes a good metal AM powder?

AM processes typically operate with 'fixed' parameters for a specific application; most currently available machines offer little opportunity for responsive control. As such, inconsistent properties in the input material will directly translate to inconsistent properties in the finished component. Poor powder quality can produce defects in the end part including pores, cracks, inclusions, residual stresses, and sub-optimal surface roughness, as well as compromising throughput. To avoid these issues, manufacturers must understand the relationship between material properties, processing performance, and end-component properties [3].

Chemistry is paramount in metal AM powders. A powder needs to be compatible with the alloy composition of the material specified. The grade must also be carefully selected to control the interstitial elements present such as oxygen or nitrogen, as well as particulate contaminants which can impact the properties of the finished part. Beyond chemistry, it is the physical characteristics of a metal powder, such as packing density and flowability, that determine AM performance. Powders that pack consistently with high density are associated with the production of consistent-quality components with fewer flaws, while good flowability enables the powder to spread evenly and smoothly across a bed to form a uniform layer with no air voids [3].

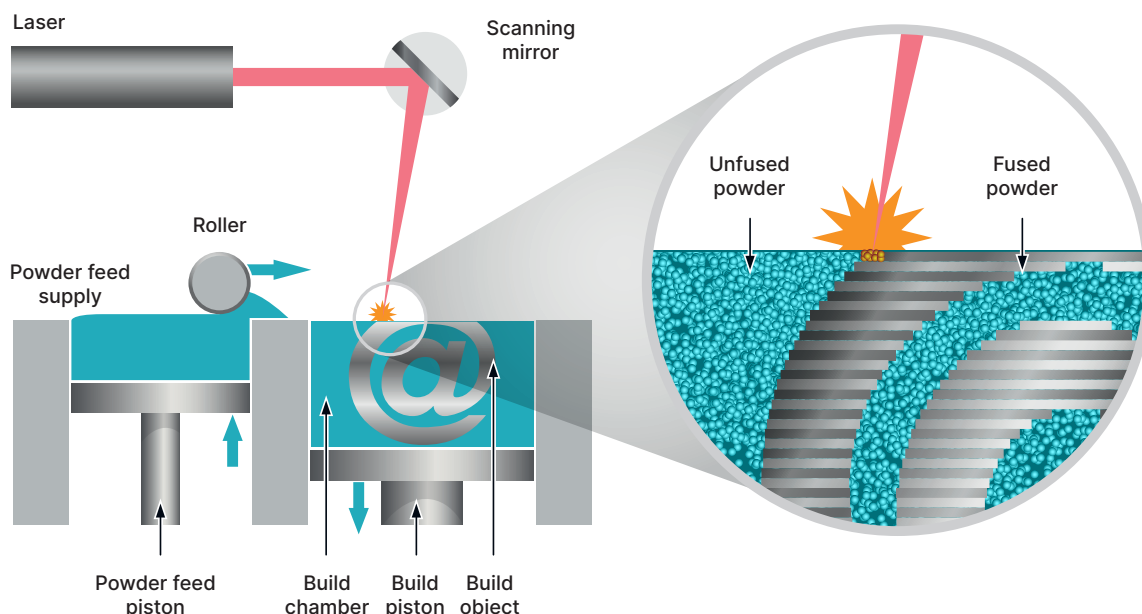


Figure 1: Powder bed AM processes such as SLM require rapid, even powder spreading and effective recycling of the excess powder.

Bulk density and flowability are directly, though not exclusively, influenced by morphological characteristics such as particle size and shape, as illustrated in Figure 2. For instance, smooth, regular-shaped particles typically flow more easily than those with a rough surface or an irregular shape [4]. This is because rougher surfaces result in increased interparticle friction, while irregularly shaped particles are more prone to mechanical interlocking; both effects decrease flowability. Similarly, spherical particles tend to pack more efficiently than those that are irregular, giving rise to higher bulk densities [5].

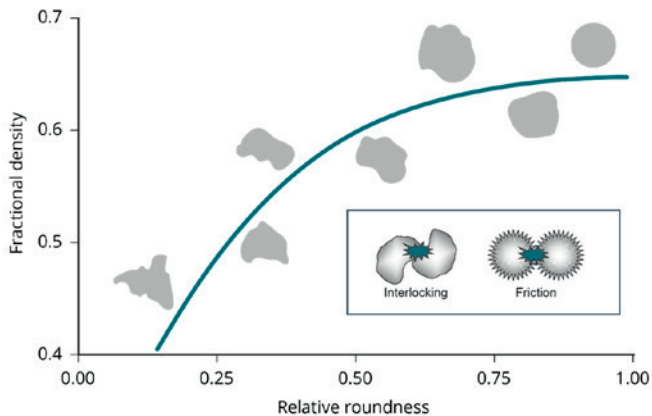


Figure 2: Smooth, regularly shaped particles tend to flow more easily than those that are irregular and/or rougher, because of reduced interparticle friction and a lower risk of mechanical interlocking.

When it comes to particle size, AM metal powders must be fine – for example, so that they can form a powder bed just tens of microns thick. However, these ‘fines’ can be problematic for health and safety, as well as flowability. Because the forces of attraction between particles increase with decreasing particle size, finer powders usually flow less freely than coarser analogs. Optimizing particle shape can help to mitigate this effect [4, 6]. Particle size and particle size distribution also influence packing, as illustrated in Figure 3. Specifically, distributions including both coarse and fine particles, with finer particles filling the interstices left by larger ones [5], enable maximum packing density.

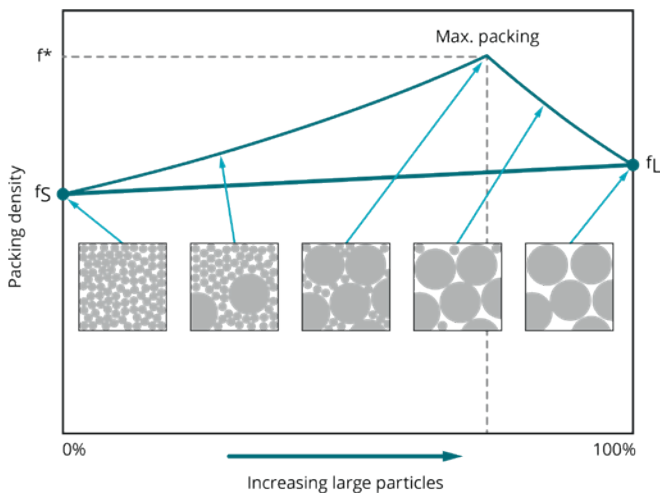


Figure 3: Packing density reaches a maximum when the particle size distribution includes both fine and coarse particles.

How does the production of metal AM powders affect their properties?

Most metal powders used in AM are produced by gas atomization. In this process, a feedstock is melted in a crucible and then ejected through a nozzle into a high-pressure gas stream (usually argon or nitrogen), breaking the molten stream into droplets. The size of particles produced by gas atomization can be controlled by varying process parameters such as gas pressure, melt properties, nozzle design, and gas-metal ratio. However, the resulting powder is not ideal for AM processes, which optimally require a narrower particle size distribution to produce a consistent powder layer of the correct thickness, as illustrated in Figure 4. Various post-atomization processes can be applied to obtain the required size fraction. These processes include ‘scalping’ to remove oversize particles, followed by either air classification or sieving. The resulting lower atomization yield is one factor that increases the cost of AM powders [3].

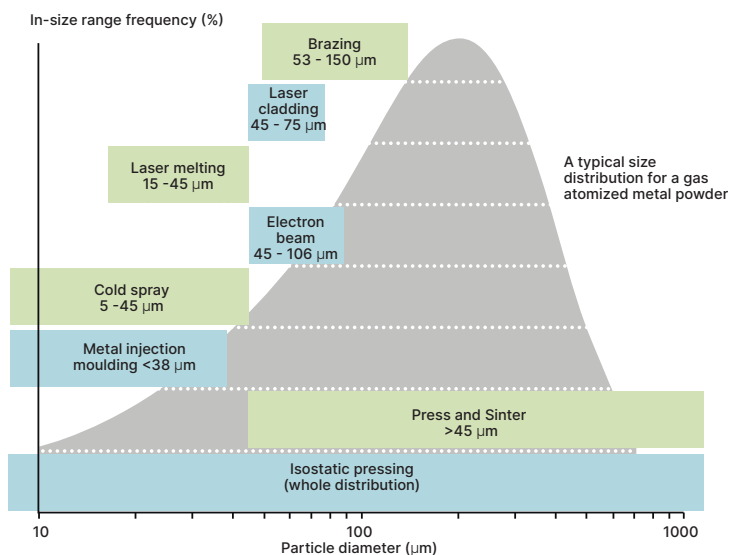


Figure 4: Typical as-atomized particle size distribution of gas-atomized powders, including required size distributions for various advanced powder metallurgy manufacturing technologies.

As stated previously, spherical particles are most suitable for powder bed AM because they have better packing and flowability. Gas-atomized particles are relatively spherical but may have unwanted features such as satellite formation, where small and larger particles fuse or agglomerate during atomization to form irregular shapes. Not only does satellite formation impair flowability and packing, but the satellite particles are so small (usually 1-10 microns) that they can also become an airborne health and safety risk if detached. More spherical particles can be produced by Plasma Atomization or the Plasma Rotating Electrode Process (PREP), but at a higher price [3].

Measuring particle size and shape

Laser diffraction

With a measurement range from 0.01 to 3,500 μm , laser diffraction is the particle sizing technology of choice for most AM applications – particularly at smaller size ranges. Laser diffraction systems determine particle size from the resulting light scattering pattern as a collimated laser beam passes through the sample, as shown in Figure 5. Large particles scatter with high intensity at narrow angles relative to the incident beam, while the signal generated by smaller particles is weaker but extends to wider angles. Laser diffraction analyzers calculate the particle size distribution of the sample from the measured angular dependence of the scattered light using an appropriate theory of light scattering, generally the Mie theory.

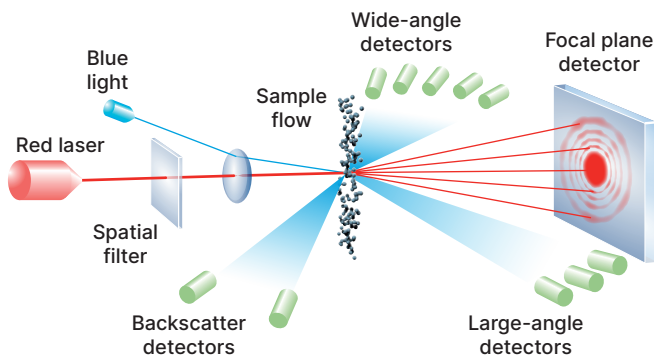


Figure 5: Illustration showing the principle behind a laser diffraction measurement, with diffracted light from dispersed particles picked up by optimally positioned detectors.

Modern laser diffraction systems such as the Mastersizer 3000 are highly automated, to the point of push-button operation, and offer high-throughput analysis with minimal manual input. Besides laboratory-based laser diffraction systems, there are also online process systems, such as Insitac, that deliver real-time particle size monitoring for automated process control. These online process systems can be used either to monitor particle size evolution during atomization, grinding, or spray drying, or at an end-user facility for automated powder handling and recycling [7].

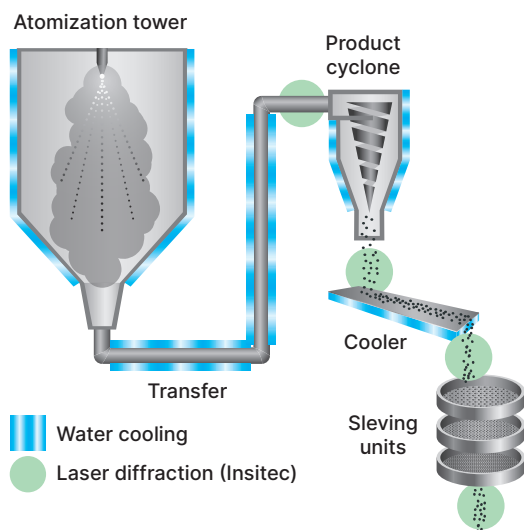


Figure 6: Schematic of a gas-atomization process for manufacturing metal powders, showing the points where laser diffraction can be employed.

Figure 7 shows measurements for four fractions of metal powders, made using both wet and dry dispersion on the Mastersizer 3000. Both wet and dry dispersion are suitable for analyzing metal powders and should give equivalent results if the dispersion procedures are optimized and sampling is comparable. For the $<150\mu\text{m}$ fraction, there is a noticeable discrepancy between the wet and dry measurements. This is likely due either to fine particles adhering to larger particles in the dry state, or a difference in sampling [8].

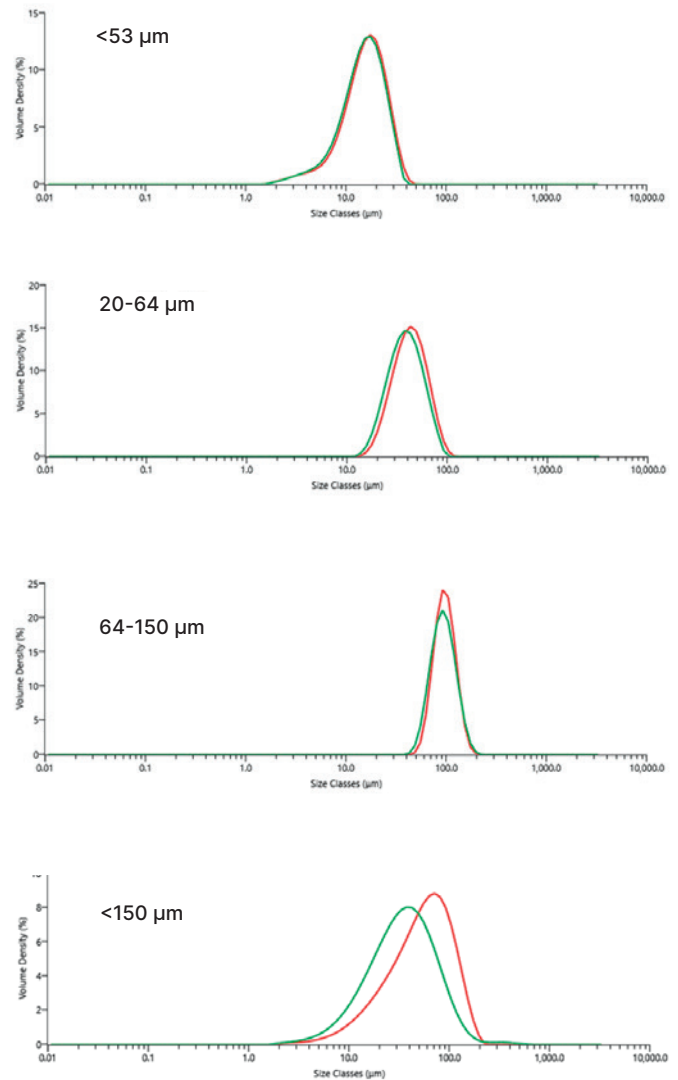


Figure 7: Comparisons of wet and dry measurements for four stainless-steel 316L powder samples. In each case, the red trace shows the dry particle size distribution (PSD) and the green trace shows the wet PSD. Each trace shows the average result over five measurements.

Table 1 below shows a comparison between dry dispersion measurements from the Mastersizer 3000 (lab analyzer) and Insitac (on-line analyzer) for four different size fractions of a stainless-steel powder. The Insitac and Mastersizer results are in good agreement for all fractions, with the Insitac results $< 2\%$ higher than the Mastersizer results [9].

Table 1: Comparison between measurements from the Mastersizer 3000 (lab analyzer) and Insitec (on-line analyzer) for four different size fractions of a stainless-steel powder.

	<25 μm	<25 μm	20-64 μm	20-64 μm	64-150 μm	64-150 μm	<150 μm	<150 μm
	MS3000	Insitec	MS3000	Insitec	MS3000	Insitec	MS3000	Insitec
d_{10}	7.46	7.11	25.3	24.9	68.9	67	16.7	16.5
d_{50}	15.9	16.5	42.3	44.2	94.7	94.2	54.7	56.7
d_{90}	27.0	27.2	68.2	70.0	131	134	117	122

Automated image analysis


When it comes to characterizing the particles used in metal AM, three main techniques are commonly used: dynamic image analysis, automated static image analysis, and scanning electron microscopy (SEM). The key distinction between these techniques is the number of individual images generated and the resolution of those images [10]:

- Dynamic image analysis yields the highest number of images, but the resolution is often too low to capture fine material or characterize surface texture.
- SEM yields the highest resolution and captures fine surface detail, but the low number of images means it can only be used as a qualitative technique.
- Static image analysis sits somewhere between SEM and dynamic image analysis, offering a resolution high enough to capture fine material and surface texture, with a statistically relevant number of images for both qualitative and quantitative analysis.

Dynamic image analysis

Dynamic imaging systems, such as the Hydro Insight, can be integrated into your laser diffraction workflow to give particle shape information at the same time as your laser diffraction measurements. Particles suspended in a flowing stream by the laser diffraction dispersion unit flow through the Hydro Insight and are then photographed by a high-resolution digital camera. The images are stored for viewing and converted to a digital format for real-time image analysis and quantification of size and/or shape.

The most significant benefits of this approach are that large numbers of particles can be measured quickly (as with laser diffraction) and that all three particle dimensions are captured, due to the random orientation of the particles. Other benefits include the ability to detect small numbers of oversize particles or contaminants that laser diffraction may miss, and to troubleshoot or optimize laser diffraction methods [11]. Additionally, the Hydro Insight also provides a built-in algorithm for correlating particle size data with sieve analysis. The algorithm measures the bounded rectangular width of irregular particles that would fit through a sieve mesh of the same size and automatically sorts the data into sieve fractions. These measurements tend to correlate better with sieve analysis than equivalent circular areas or spherical diameters reported by laser diffraction, especially when dealing with non-spherical particles [12].

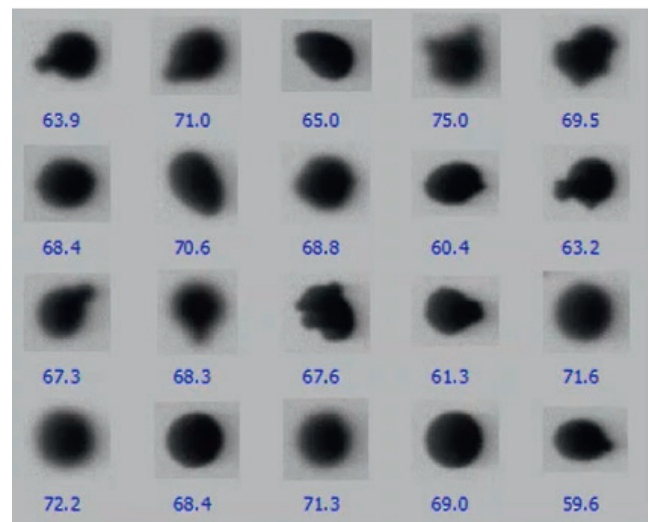

Aperture size (μm)	ECA diameter %	BR width %	Differences
<20	0.02	0.543	-0.523
20	2.017	6.378	-4.361
25	16.818	24.692	-7.874
32	22.987	26.327	-3.34
38	23.495	15.666	7.829
45	9.817	2.496	7.321
53	3.913	3.071	0.842
63	3.265	3.497	-0.232
73	4.254	4.585	-0.331
90	2.257	1.825	0.432
103	2.747	2.450	0.297
125	4.217	3.688	0.529
150	2.280	2.772	-0.492
180	1.913	2.011	-0.098

Figure 8: Thumbnail images of metal powders, captured by Hydro Insight (top), and sieve correlations based on equivalent circular area (ECA) diameter and bounded rectangle (BR) width (bottom).

Static image analysis

Suitable for particles from approximately $0.5\ \mu\text{m}$ to $> 1\ \text{mm}$, static imaging systems such as Morphologi 4 capture individual images of tens of thousands of particles. From these images, multiple size and shape parameters can be calculated. Unlike in dynamic imaging, the sample is dispersed on a substrate and remains effectively static. The substrate is then moved relative to the microscope optics, situated above the sample, using a motorized stage. The sample is illuminated from either below the stage (diascopic) or above it (episcopic), depending on the image requirements. The captured images are processed for size and shape analysis. Figure 9 illustrates this workflow.

Because the distance between the microscope and substrate is well controlled, all the particles remain in focus during the measurement, so the resulting images are of higher quality than with a dynamic approach. Static imaging also reliably captures the two largest particle dimensions of irregular particles, since they take a preferred orientation on the substrate. Furthermore, because the position of each particle is recorded, particles of interest can be revisited with higher magnification for a more detailed study.

The Morphologi 4 provides 20 different shape parameters, including circularity, elongation, convexity, and solidity. It can also support customized classifications to examine features such as satelliting. Figure 10 shows several metal powder images taken with the Morphologi 4. The instrument has automatically classified and grouped these images based on their shape, taking into account parameters such as sphericity, elongation, and the presence of satellites [13].

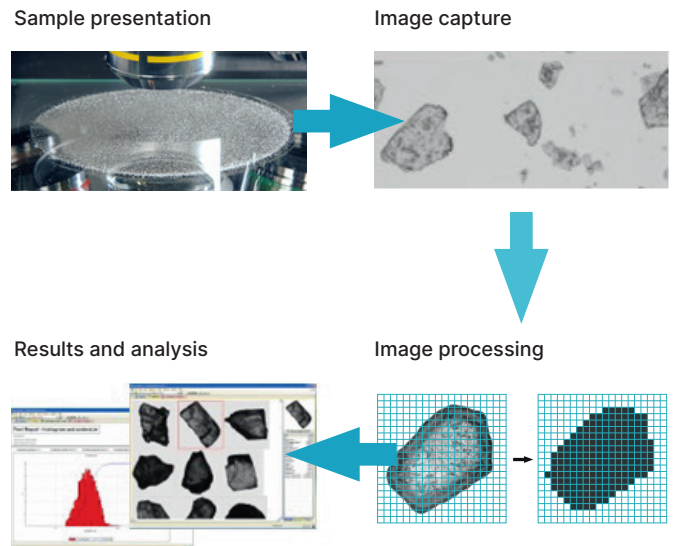


Figure 9: Illustration showing the general workflow for an automated static imaging measurement.

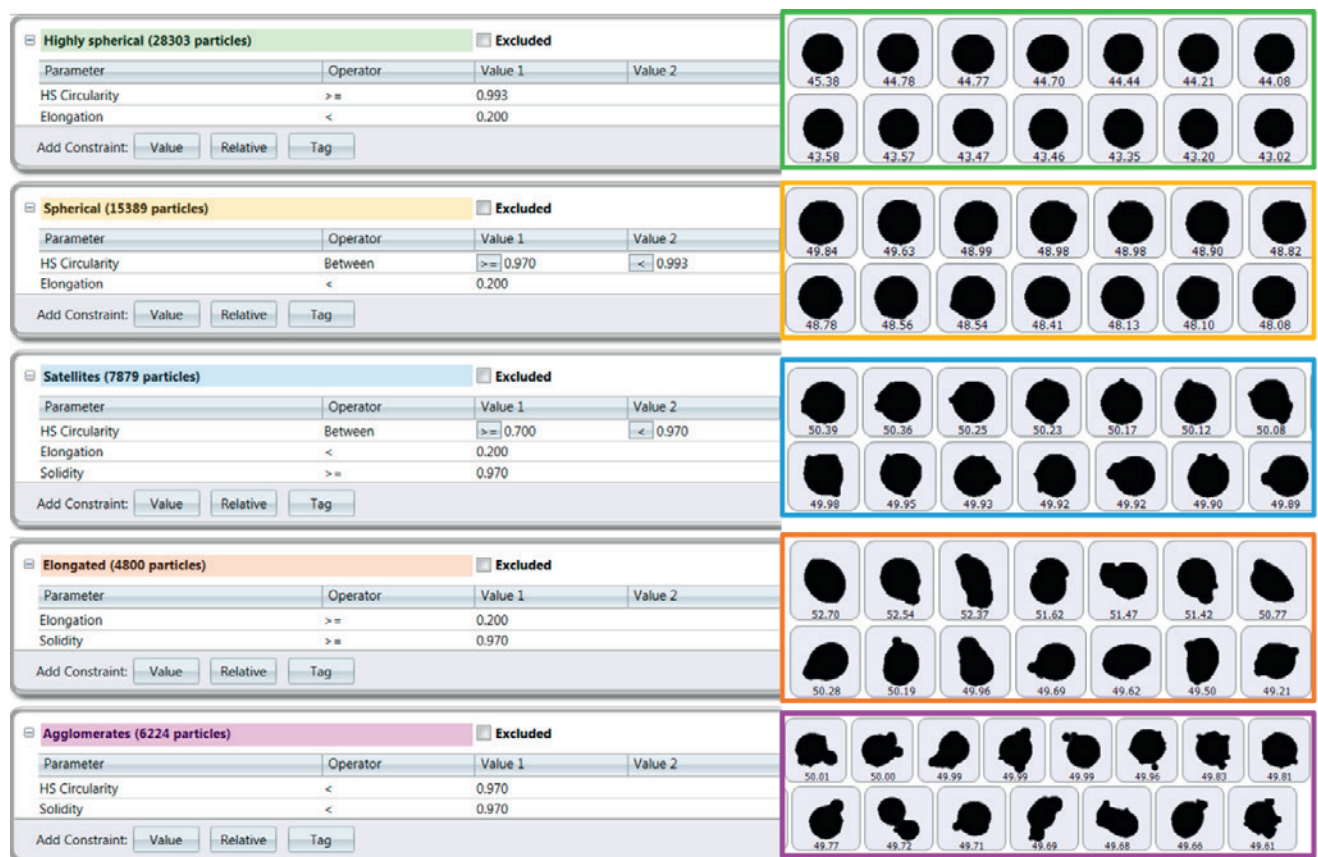


Figure 10: Particle classifications for an AM metal powder with corresponding particle images.

Conclusions

The size and shape of metal powder particles affect powder bed packing and flowability. In turn, these features impact the build quality and final properties of components manufactured using AM. As such, understanding and optimizing particle size and shape is critical to the success of powder-bed AM. Laser diffraction and automated image analysis are complementary tools that can be used to characterize and optimize metal powders for a range of powder-bed AM processes.

- [1] Wohlers Report 2020. 3D Printing and Additive Manufacturing State of the Industry. Annual Worldwide Progress Report
- [2] SmarTech Publishing 'The Top Three additive Manufacturing Predictions for 3D printing in 2017'
- [3] The importance of powder quality in powder bed Additive Manufacturing processes: Malvern Panalytical webinar available for viewing
- [4] D.F. Heaney, Handbook of metal injection molding, Woodhead Publishing, 2012
- [5] J.P. Bennett & J.D. Smith, Fundamentals of Refractory Technology (Ceramic Transaction Series), Volume 25, 2001 (American Chemical Society)
- [6] C.N. Davies, Aerosol Science, Academic Press, London and New York, 1966
- [7] J. DeNigris, "Taking control of metal powder properties: Exploring the benefits of real-time particle sizing," Met. Powder Rep., 73, No. 4, 202–207 (2018)
- [8] Determining the particle size distribution of metal powders using wet and dry dispersion on the Mastersizer 3000: Malvern Panalytical Application Note available for download
- [9] M. Tulley, S. Hall, U. M. Attia, J. Dawes, J. Ashby, and G. Thornton, Feasibility Assessment of Using In-Process Measurement Analysers for Metal Powders, Euro PM
- [10] 8 reasons why it's time to upgrade to automated imaging: A Malvern Panalytical Application Note available for download
- [11] Five compelling reasons to add the Hydro Insight dynamic imaging accessory to your Mastersizer: Malvern Panalytical White Paper available for download
- [12] A complete solution for particle size analysis of metal powders - A combination of laser diffraction and imaging: Malvern Panalytical Application Note available for download
- [13] Characterising the particle size and shape of metal powders for Additive Layer Manufacturing: Malvern Panalytical Application Note available for download

Why choose us?

**When you make the invisible visible,
the impossible is possible.**

Our analytical systems and services help our customers to create a better world. Through chemical, physical and structural analysis of materials, they improve everything from the energies that power us and the materials we build with, to the medicines that cure us and the foods we enjoy.

We partner with many of the world's biggest companies, universities and research organizations. They value us not only for the power of our solutions, but also for the depth of our expertise, collaboration and integrity.

With over 2200 employees, we serve the world, and we are part of Spectris plc, the world-leading precision measurements group.

Malvern Panalytical. We're BIG on small™

Service & Support

Malvern Panalytical provides the global training, service and support you need to continuously drive your analytical processes at the highest level. We help you increase the return on your investment with us, and ensure that as your laboratory and analytical needs grow, we are there to support you.

Our worldwide team of specialists adds value to your business processes by ensuring applications expertise, rapid response and maximum instrument uptime.

- Local and remote support
- Full and flexible range of support agreements
- Compliance and validation support
- Onsite or classroom-based training courses
- e-Learning training courses and web seminars
- Sample and application consultancy



 **Prager
Elektronik**

Traunstraße 21, A-2120 Wolkersdorf
T: +43 2245 6725 F: +43 2245 559633
office@prager-elektronik.at
www.prager-elektronik.at

Malvern Panalytical

Grovewood Road, Malvern,
Worcestershire, WR14 1XZ,
United Kingdom

Tel. +44 1684 892456
Fax. +44 1684 892789

Lelyweg 1,
7602 EA Almelo,
The Netherlands

Tel. +31 546 534 444
Fax. +31 546 534 598